# Q-Sun Operator Training

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[View Recorded Presentation](#)

Presented By

Chanatorn Chantanukul  
Technical Sales of H.J.Unkel (Thai) Limited

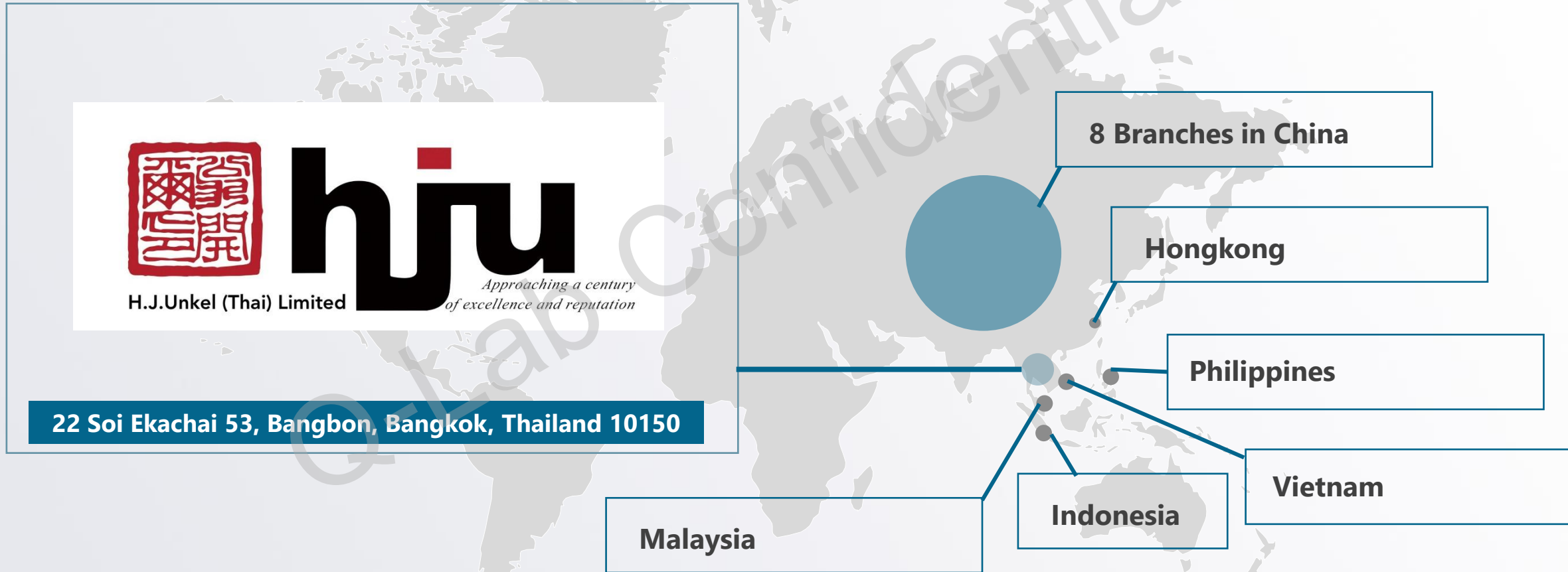
H.J.Unkel (Thai) Limited

*Approaching a century  
of excellence and reputation*

# H.J.UNKEL GROUP

Serve Quality technical and chemical products. throughout China and Southeast Asia

**hju**  
Approaching a century  
of excellence and reputation

H.J.Unkel (Thai) Limited

# 47 Years Experience

Serve Quality technical

1976

"Helmut & Associated Limited" was established

at Ratchadamri,  
Bangkok

1990

"Helmut & Associated Limited"

changed name to

"H.J. Unkel (Thailand) Limited"

1993

"H.J. Unkel (Thailand) Limited"

changed name to

"H.J. Unkel (Technic) Limited "

2007

"H.J. Unkel (Technic) Limited"

changed name to

"H.J.Unkel (Thai) Limited"

2020

"H.J. Unkel (Thai) Limited" have finally gotten ISO 9001:2015 Certificate issued by TISTR.

# 35 Years cooperation With Q-Lab

Q-Lab Corporation is a global provider of material durability testing products since 1956. We design and manufacture standard test substrates as well as weathering, light stability, and corrosion testers.

# 35 Years cooperation

## With Q-Lab

▲ Q-Lab Representative Meeting in USA

▲ Q-Lab Sales Training in Shanghai

# We Have The Best Team

## HEAD OF SALES - EQUIPMENT

### PROFILE

Hold a Bachelor of Business Administration Program in Business Computer

### EXPERIENCE

15 years in sales at H.J.Unkel

**CHANATORN C.**

## HEAD OF TECHNICAL SERVICE

### PROFILE

Hold a Technical Diploma in Industrial Technology; specialized in Electrical Power & Installation

### EXPERIENCE

17 years in technical field  
12 years with H.J.Unkel

**SUCHART S.**

## CUSTOMER TECHNICAL SUPPORT

### PROFILE

Hold a Bachelor of Science Program in Production Technology Management; specialized in PLC

### EXPERIENCE

13 years in technical field at H.J.Unkel

**ANUSORN S.**

# Q-Lab Corporation

- Founded in 1956
- Specialize in material durability testing equipment and services

**Westlake, Ohio  
Headquarters &  
Instrument Division**

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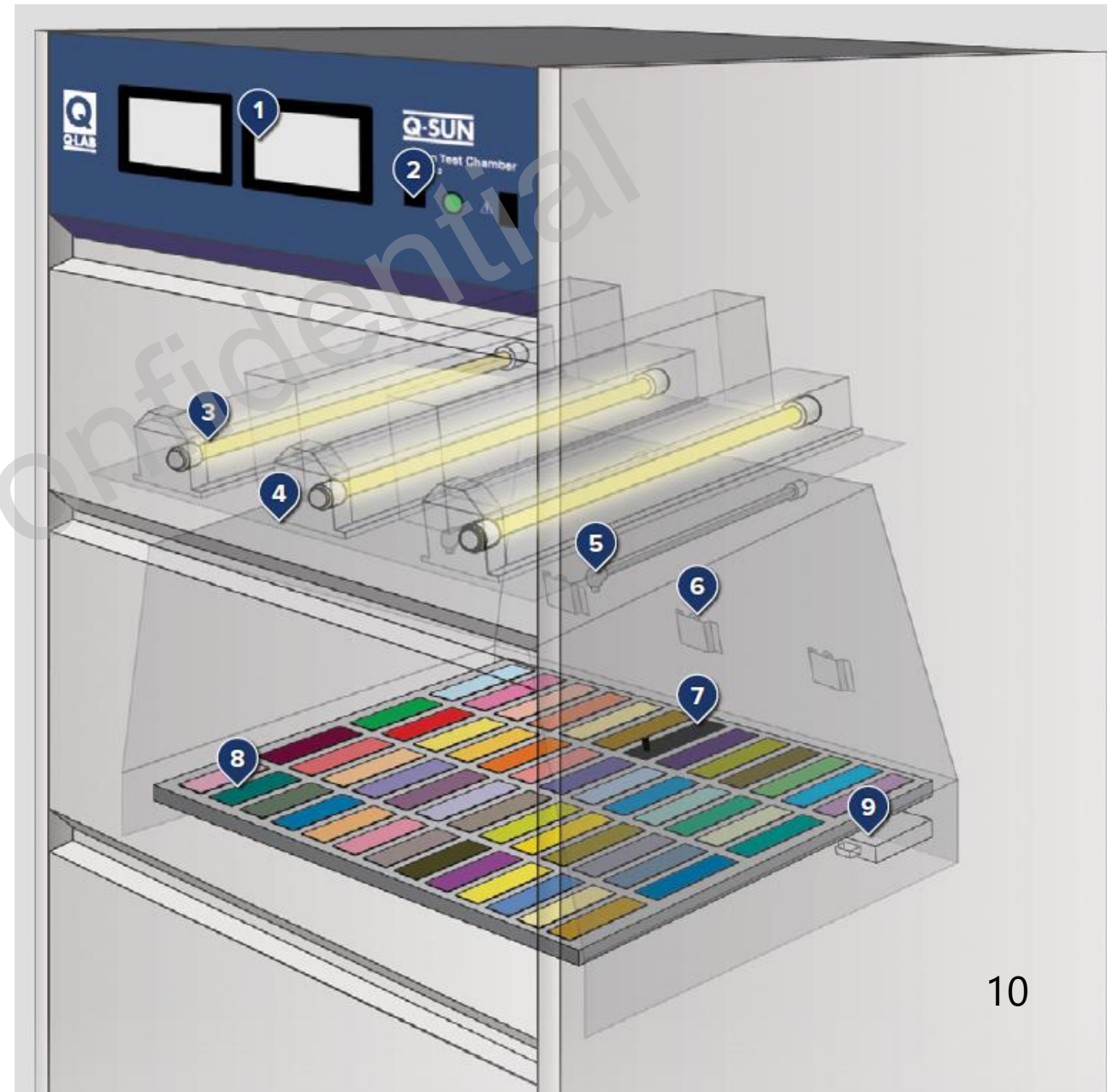
**Xe-2**


# Flat Array

- 1) User interface
- 2) USB port for data transfer
- 3) Xenon lamps with irradiance control
- 4) Optical filters
- 5) Water spray
- 6) Onboard irradiance sensors
- 7) Black Panel Temp sensor
- 8) Specimen holders
- 9) Relative Humidity/CAT sensor

# Rotating Rack

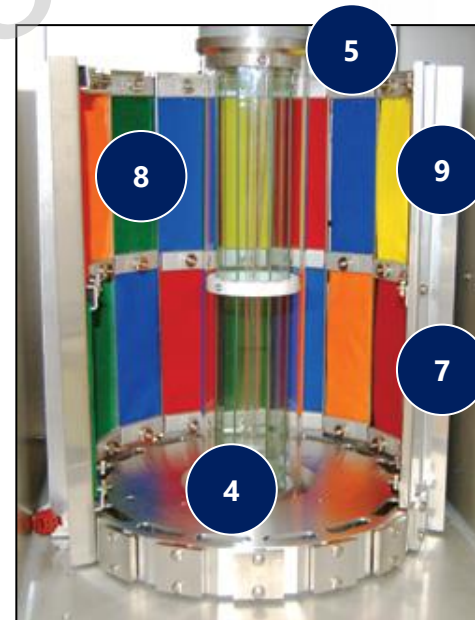- 1) User interface
- 2) USB port for data transfer
- 3) Xenon lamps with irradiance control
- 4) Optical filters
- 5) Water spray
- 6) Onboard irradiance sensors
- 7) Black Panel Temp sensor
- 8) Specimen holders
- 9) Relative Humidity/CAT sensor

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Model/ Configuration		Tester Features					
		Gen 4 + High Irradiance (-E) <sup>3</sup>	Humid Control (-H)	Chamber Air Chiller (-C)	Water Spray (-S)	Back Spray (-BS) <sup>6</sup>	Dual Spray (-DS) <sup>6</sup>
Xe-1	Xe-1-BE	●					
	Xe-1-BCE	●		●			
	Xe-1-SE	●			●		
	Xe-1-SCE	●		●	●		
	Xe-1-WE	●			● <sup>4</sup>		

Xe-2	Xe-2-HE	●	●				
	Xe-2-HSE	●	●		●		
	Xe-2-HBSE	●	●		●	●	

Xe-3	Xe-3-HE	●	●				
	Xe-3-HCE	●	●	●			
	Xe-3-HSE	●	●		●		
	Xe-3-HSCE	●	●	●	●		
	Xe-3-HBSE	●	●		●	●	
	Xe-3-HBSCE	●	●	●	●	●	
	Xe-3-HDSE	●	●		●		●
	Xe-3-HDSCE	●	●	●	●		●
	Xe-3-HDSBSE	●	●		●	●	●
	Xe-3-HDSBSCE	●	●	●	●	●	●

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
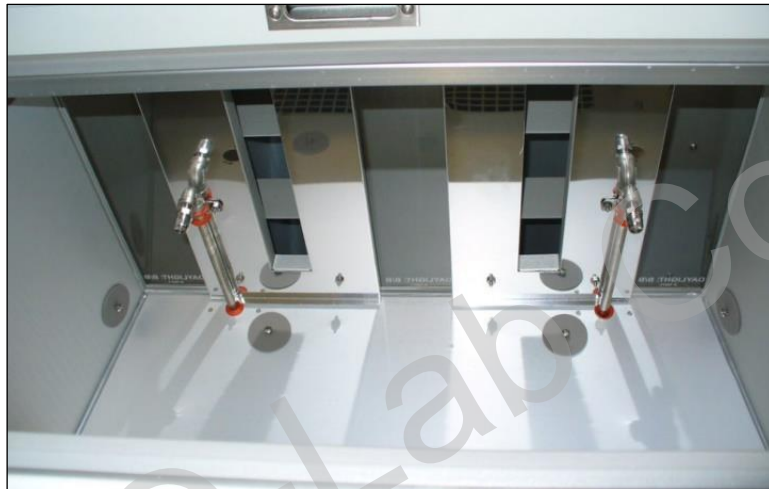
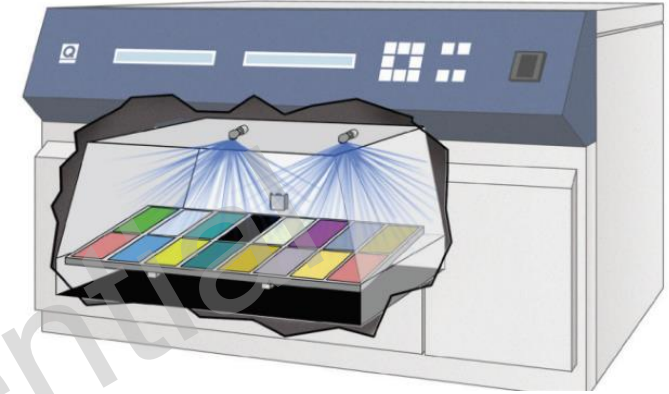
**Xe-1**

### Features

- Pulse Rate Control
- Automatic Fault Detector

### Two nozzles

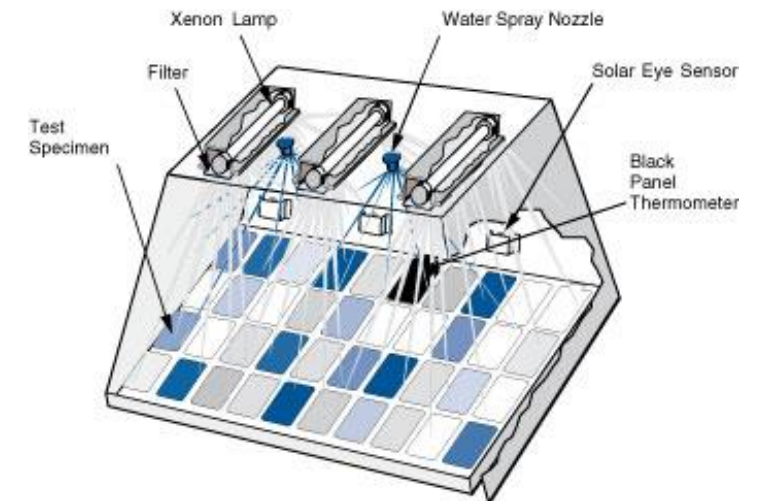- Mist
- Thermal Shock
- Erosion

**Xe-3**

# WATER SPRAY

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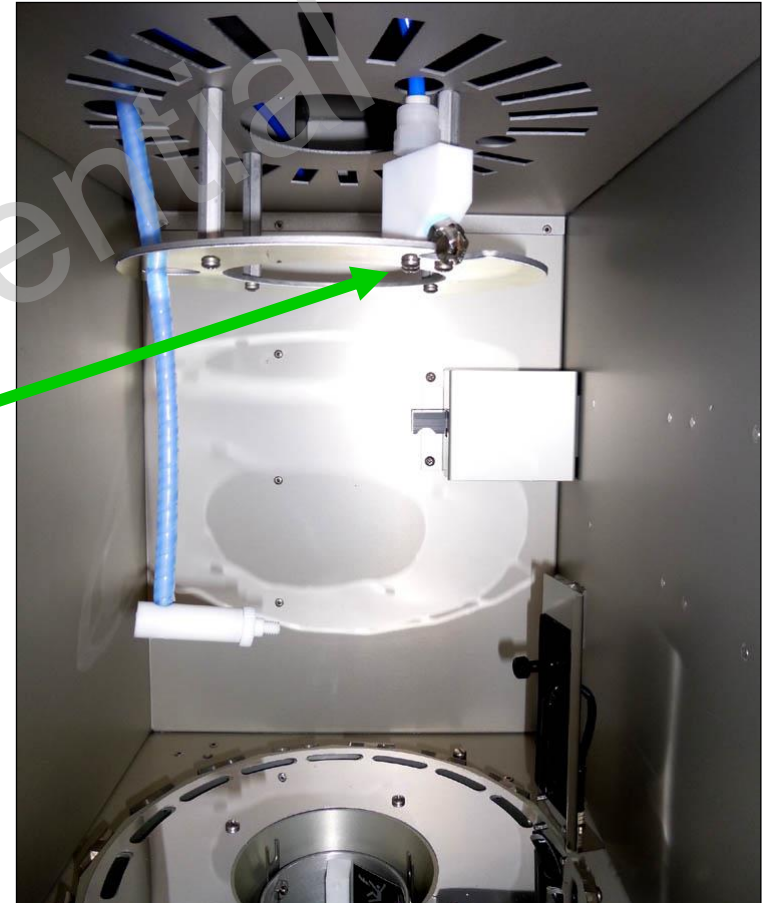
### Features

- Pulse Rate Control
- Automatic Fault Detector

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- Feedback Loop System
- RH/CAT Sensor
- Main Controller
- Humidity Generator
  - Nebulizer in Xe-2
  - Boiler in Xe-3

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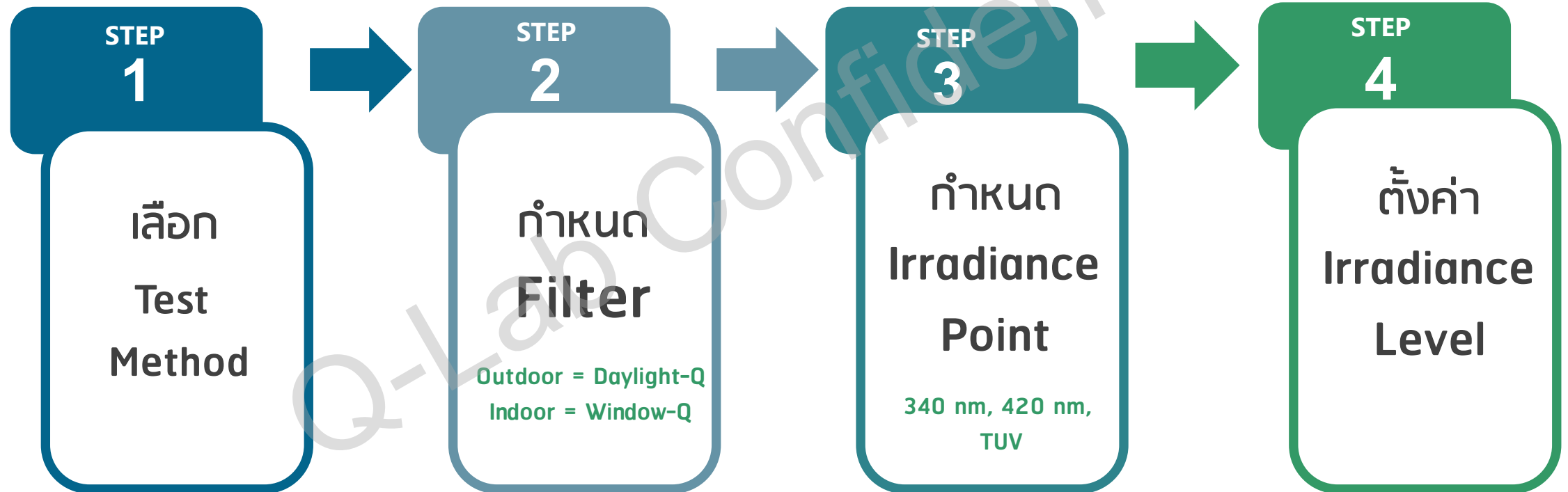
while the Xe-2 filter lantern consists of an outer borosilicate or quartz cylinder and 14 inner filters.


	Xe-1 & Xe-3 Irradiance Values Typical (& Maximum) <sup>A,B,C</sup>			Xe-2 Irradiance Values Typical (& Maximum) <sup>A,B,C</sup>				
	W/m <sup>2</sup> /nm @ 340 nm	W/m <sup>2</sup> /nm @ 420 nm	W/m <sup>2</sup> @TUV (300-400 nm)	W/m <sup>2</sup> /nm @ 340 nm	W/m <sup>2</sup> /nm @ 420 nm	W/m <sup>2</sup> @TUV (300-400 nm)		
Daylight-F	0.80 (1.30)	1.50 (2.40)	75 (125)	0.80 (0.95)	1.50 (1.70)	75 (85)		
Daylight Q	0.68 (1.10)			0.68 (0.80)				
Extended UV (-Q/B, -Quartz <sup>D</sup> )				0.51 (0.61) <sup>E</sup>				
Daylight-B/B				70 (108)			0.55 (0.65)	
Window (-Q, -B/SL)	0.55 (0.85)			42 (68)			-	42 (62)
Window (-SF5, -IR, -B04 <sup>F</sup> )	-							

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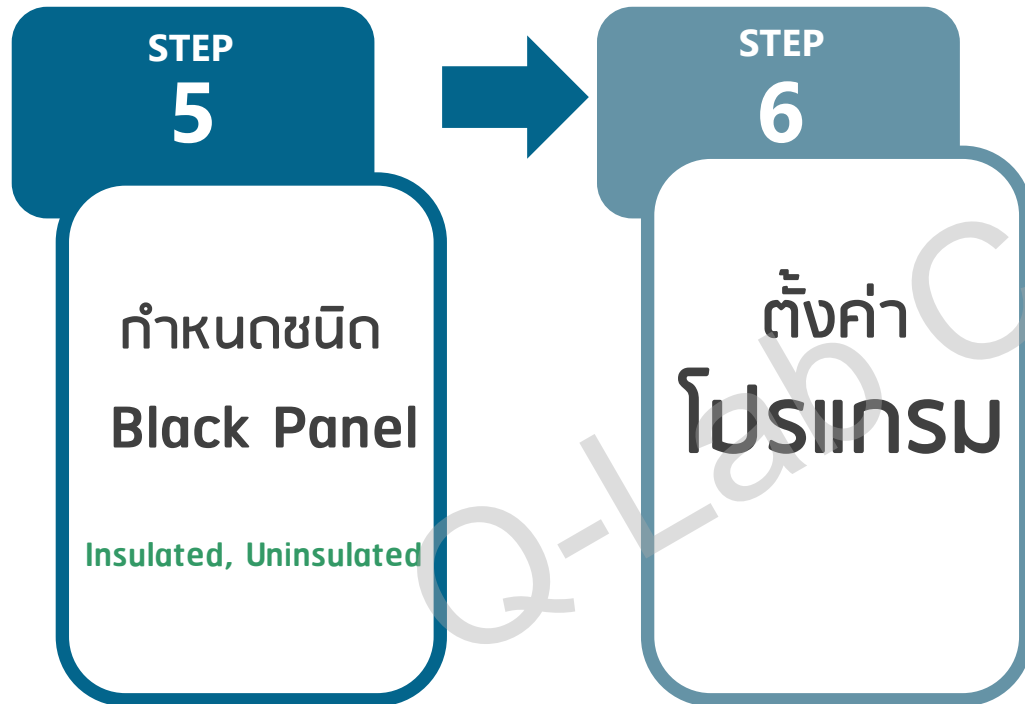## Programming

- Light
- Dark
- Light + Spray
- Dark + Spray
- Dark + Spray Front and Back
- Dual Spray
- Light + Dual Spray
- Light + Immersion
- Dark + Immersion

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**Dual, Full-Color Touchscreen Displays**

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Status Icon      State      Cycle ID and Test Name

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
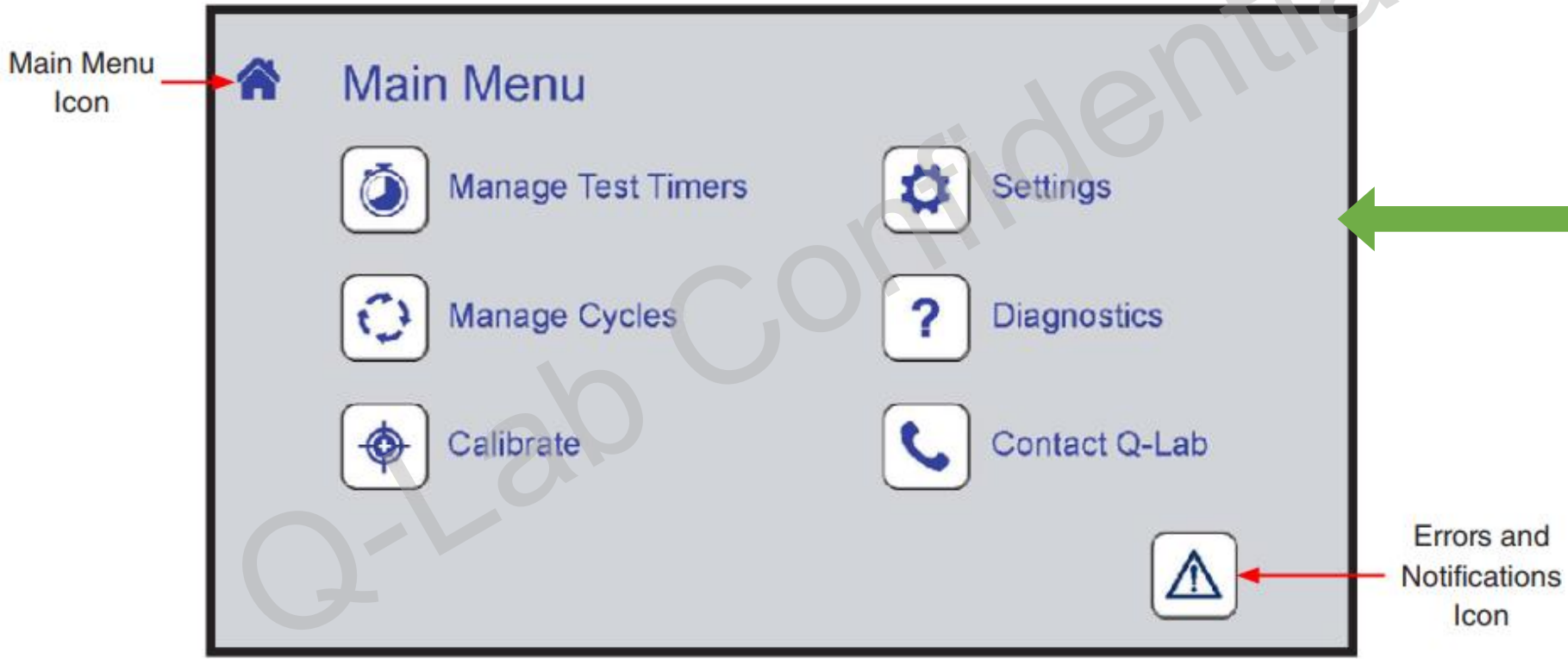
# Programing

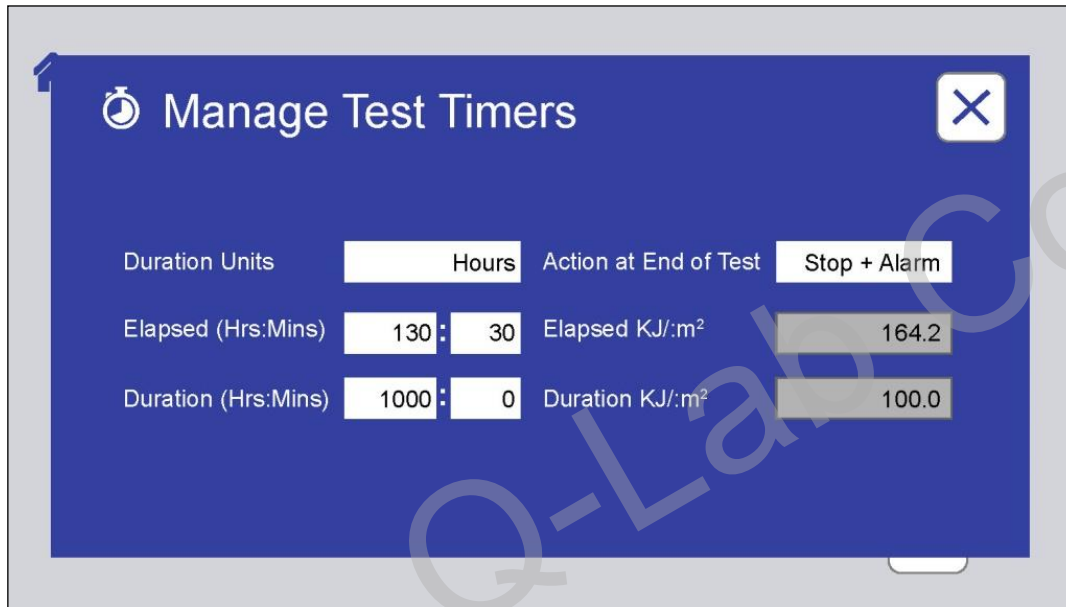
**Manage Test Timers**

Duration Units:  Action at End of Test:

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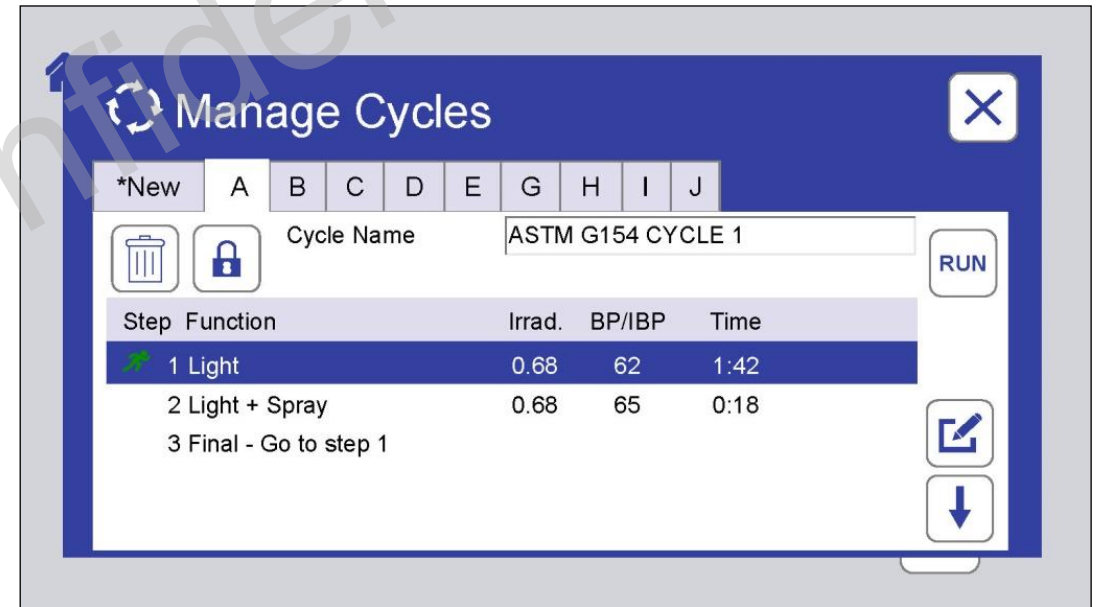- Xe-1: 17 (51 × 102 mm)
- Xe-2: 31 (45 × 132 mm)
- Xe-3: 55 (51 × 102 mm)

### Type of holder

- Open-Backed (thick, rigid specimens)
- Solid-Backed (flexible specimens)
- Three-dimensional

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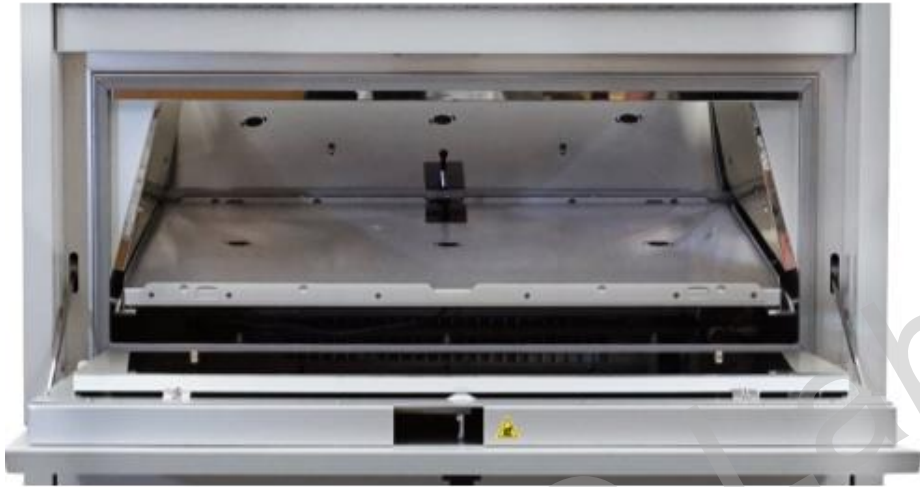
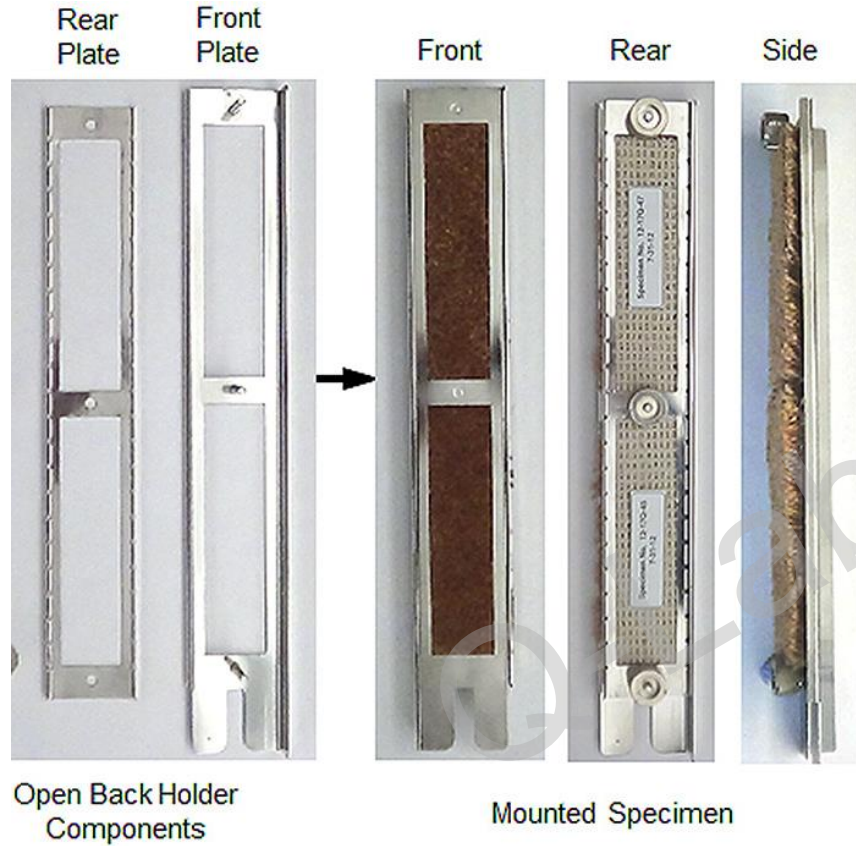
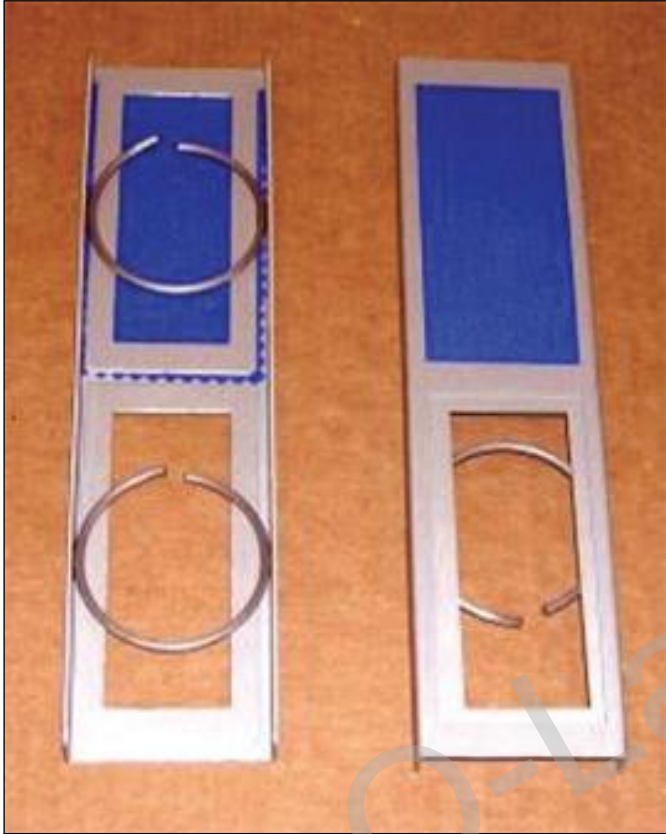
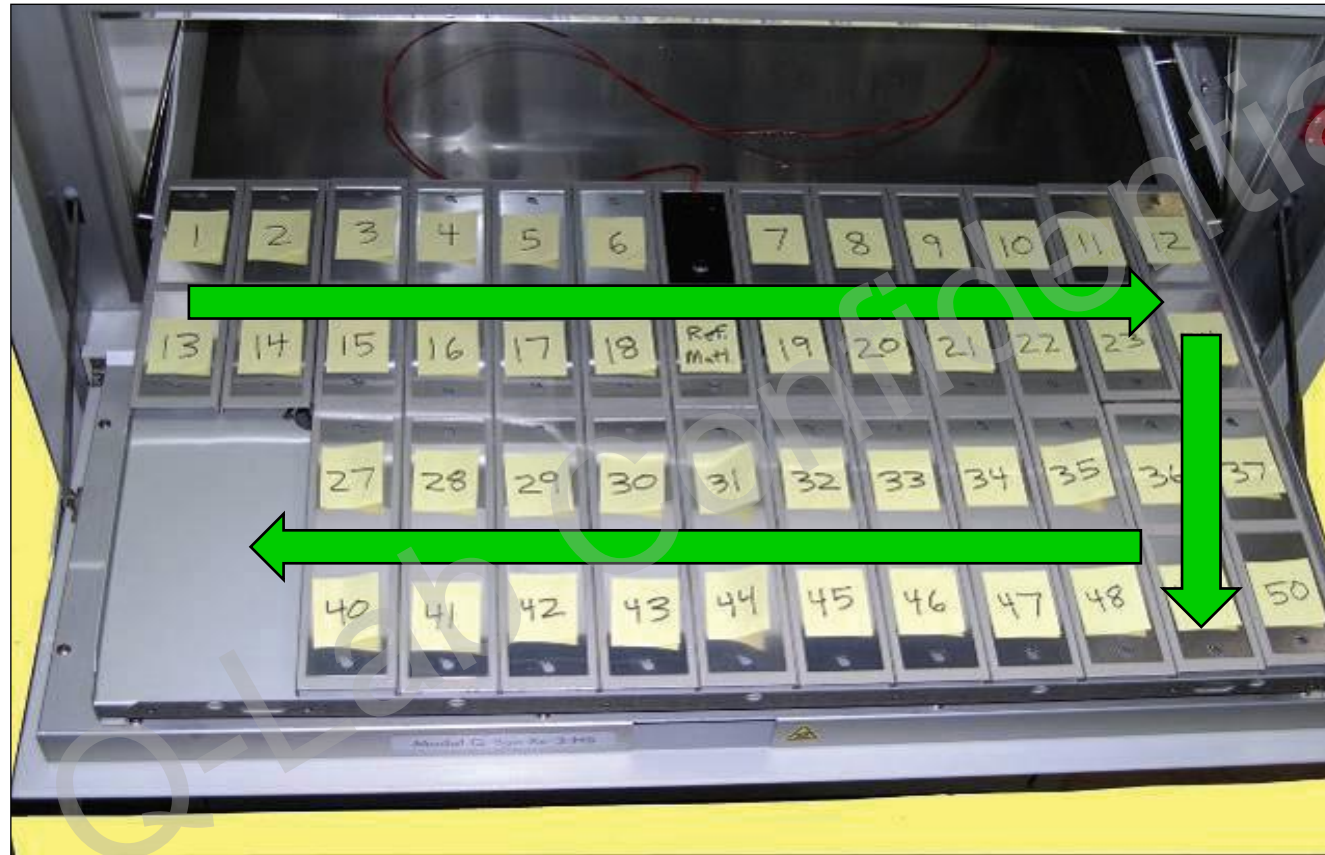
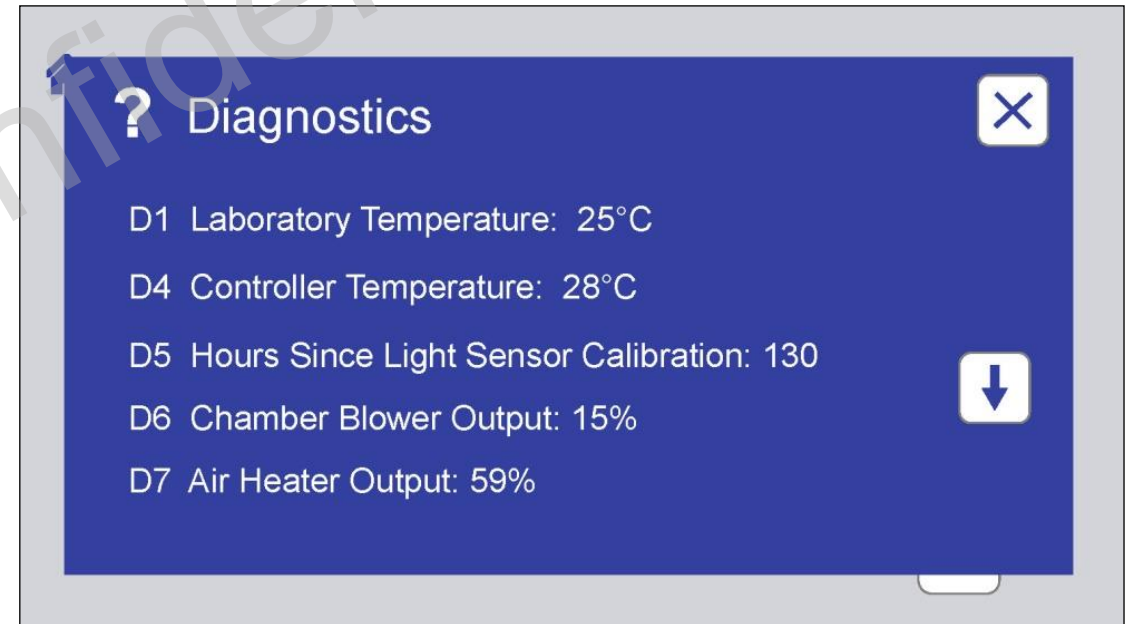
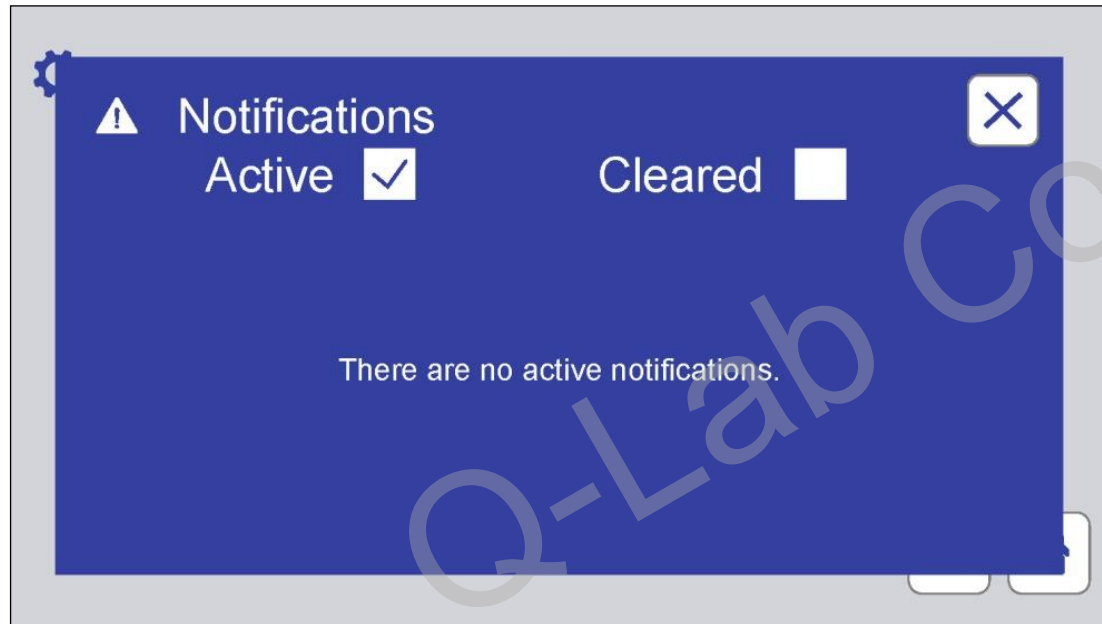
## Irradiance

UC20/340

Description	Name	Color Code	Used For
Irradiance Smart Sensor	UC20/340	Green	Calibration Radiometer for 340nm on board sensor
	UC20/420	Red	Calibration Radiometer for 420nm on board sensor
	UC20/TUV	Grey	Calibration Radiometer for TUV on board sensor
	UC20/LUX	Grey	Calibration Radiometer for LUX on board sensor

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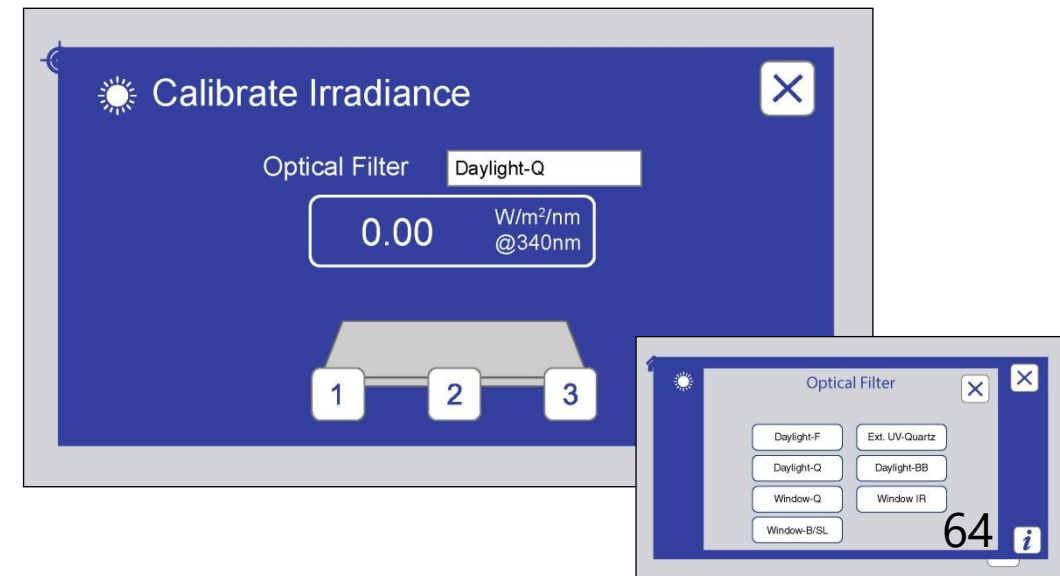
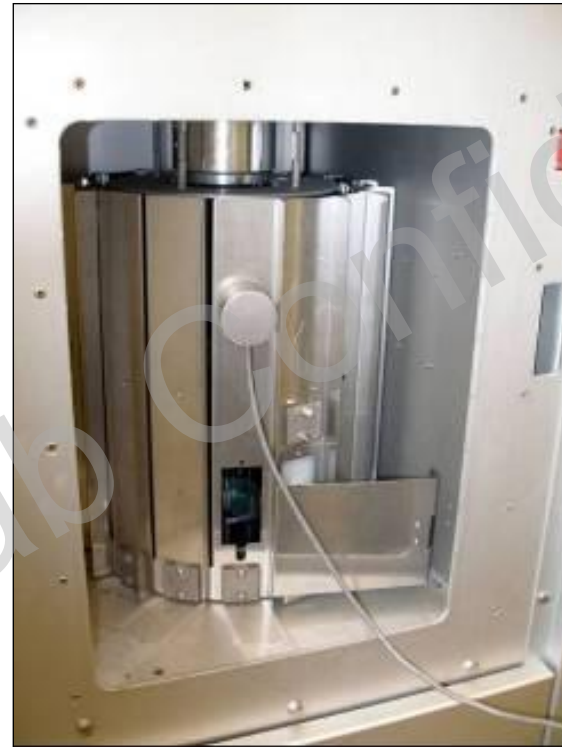
# Smart Sensor

## Temperature

Description	Name	Color Code	Used For
Temperature Smart Sensor	UC202/BP		Calibration of onboard Black Panel sensor
	UC202/IBP		Calibration of onboard Insulated Black Panel sensor

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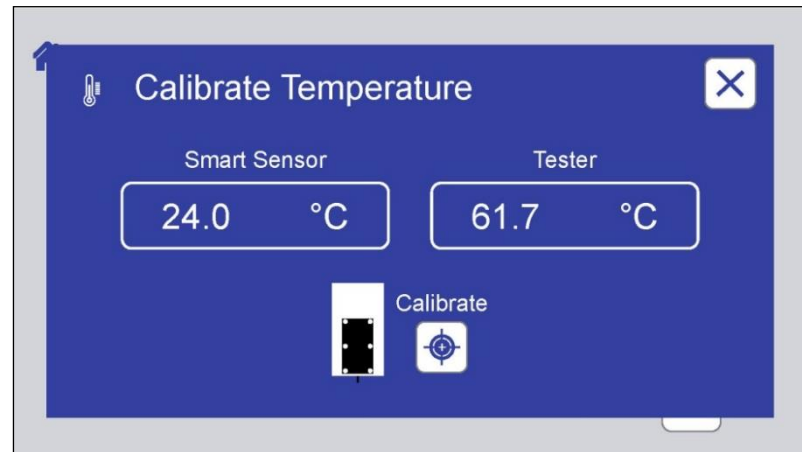
# Black Panel Temp Sensor Calibration

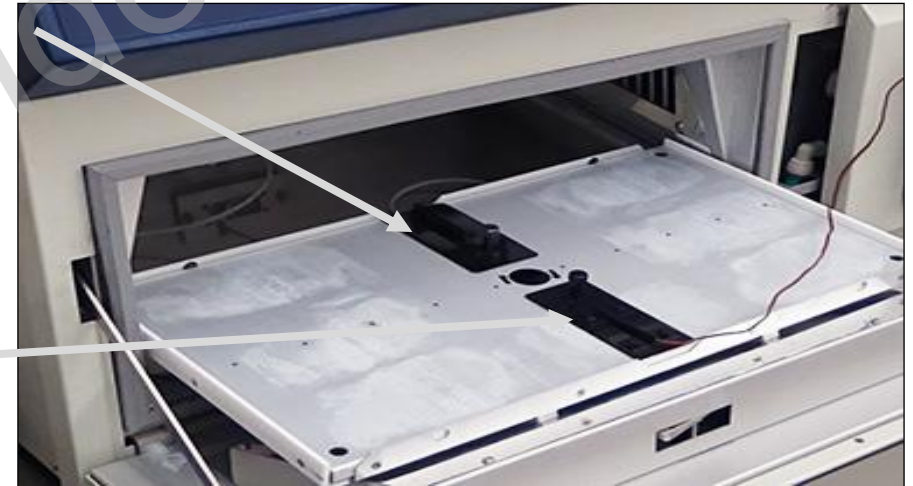
Temperature

**Onboard Sensor in Panel Holder**

Onboard Black Panel  
Calibration Black Panel

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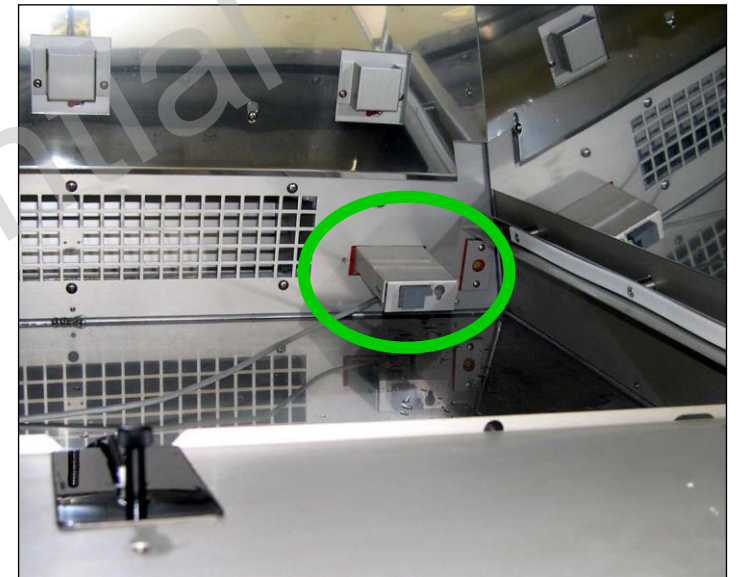
Q-SUN tester	Irradiance	Warranted Lamp Life (hrs)
Legacy	Typical	1500
Q-SUN "E" Model	Typical	3000
Q-SUN "E" Model	Maximum	1000

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